## **FOSS**

# Kjeltec™ 8200 distillation unit

### Automated distillation for Kjeldahl analysis in food and feed







Kjeltec<sup>™</sup> 8200 is a fully automated distillation unit designed primarily for Kjeldahl analysis, but widely used for many different distillation chemistries.

#### A versatile solution that reduces operational costs

A self-adjusting cooling water system saves water and reduces cost, while long-life components ensure minimal downtime. Corrosion-free materials and alkali resistant plastic splash head/tube draining vessel secure long instrument lifetime.

#### Innovative features and built in safety

Ensure safe operations with patented SAfE feature for safe distillation of samples that form salt cakes and bellow pumps for accurate reagent dispensing. A variable output steam generator broadens the application area to other volatile components.

#### Secured for your future needs

The Kjeltec<sup>™</sup> 8200 can be connected to an optional external titrator. As your business grows the Kjeltec can be upgraded to a fully automated Kjeltec 8400 analyser with a colorimetric titration system and an optional 20 or 60 place sampler.

#### Sample type

Raw materials and finished products in food, feed and agriculture

#### **Parameters**

Nitrogen, protein

#### **Technology**

Fully automated distillation for Kjeldahl and other nitrogen based chemical analysis.

#### Approvals

Officially approved (AOAC, EPA, DIN, ISO) for simplified validation

## Specifications

Performance data at 230 V	
Analysis time	3.5 minutes at 30 mg N (6.5 minutes at 200 mg N)
Distillation capacity	~ 40 ml/min at 230 V (at 100 % steam addition)
Measuring range	0.1 – 210 mg N
Reproducibility	1% RSD (including the digestion step)
Recovery	> 99.5% at nitrogen levels between 1 – 200 mg N
Tube draining	200 ml in less than 10 seconds
Reagent pump volume	0 - 150 ml in steps of 10 ml
Delay	0 – 1800 seconds
SAfE (Patented)	0 – 15 seconds
Variable steam addition	30-100 % in steps of 10 %

Installation requirements:	
Power supply	200 – 240 V 50 – 60 Hz
Tolerable voltage variation	± 10 %
Power consumption	2 200 W
Water consumption (during distillation only)	3 I/min at water temp of 25°C* 1 I/min at water temp of 15°C *Circulating Cooler recommended at >25°C
Ambient temperature	5 - 40 °C
Ambient humidity	max. 80% relative
Drain	For cooling water only
Weight net	35 kg
Dimensions (w×d×h) cm	48 × 58 × 69

#### Accessories

- Titration unit complete, 50 ml
- Basket for 20 receiver flasks (250 ml)
- Receiver flasks, 250 ml (set of 10)
- Kit for distillation by-pass
- Cable for connection of external titrator
- Kit for uprade of Kjeltec 8200 Distillation Unit to Kjeltec 8400 Analyser Unit